



# Pairing your computer to the ErgoBot: **Android**


1. Turn on the ErgoBot using the switch on its bottom.

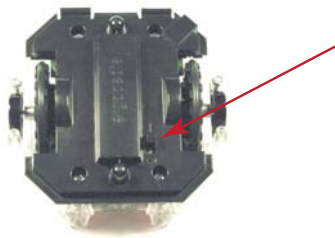


2. On your Android OS device, swipe your finger down in the upper-right corner to see the menu.

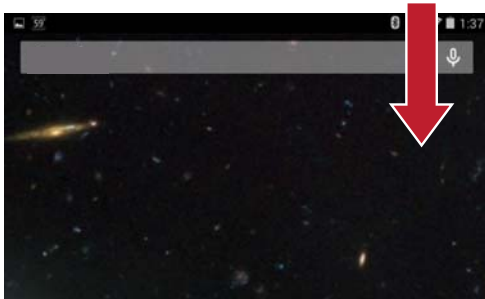


3. Select “Bluetooth” from the menu. (Either a single press or a long press, depending on your device.)

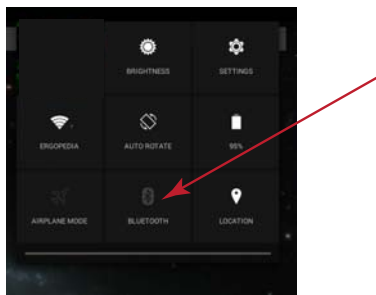


4. Turn Bluetooth On.

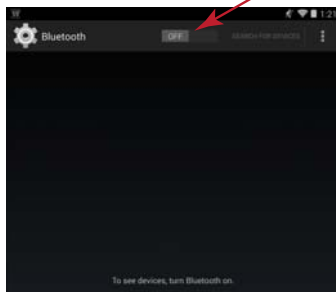


5. Select “SEARCH FOR DEVICES” to add the new ErgoBot as a Bluetooth device to your mobile device.

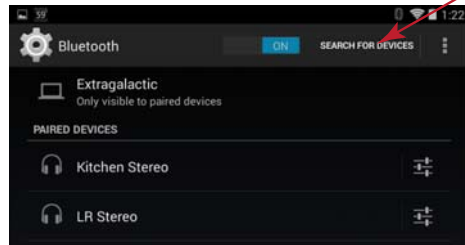

6. Select the ErgoBot device to which you want to pair and connect.

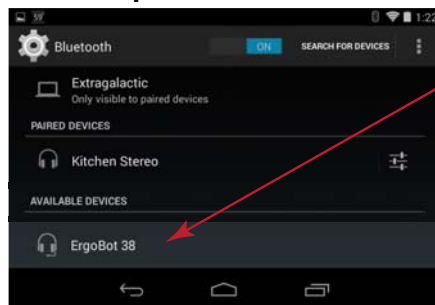

7. This shows a successful connection.

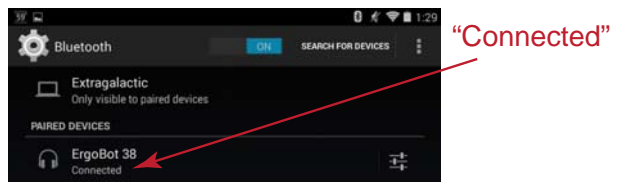

8. Sometimes you may have to turn the ErgoBot power switch on and off to refresh the Bluetooth connection.

You can now run the interactive elements in the *Essential Physics* e-Book with your paired ErgoBot!