



Pairing your computer to the ErgoBot: **Chromebook**


1. Turn on the ErgoBot using the switch on its bottom.

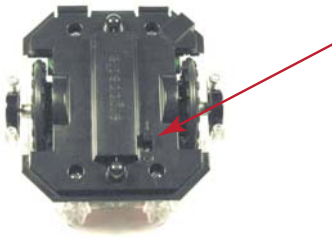


2. On your Chromebook, select the system tray in the lower left corner.

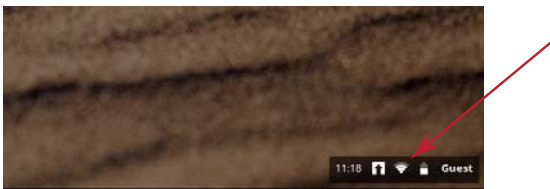

3. Select "Bluetooth" from the menu.

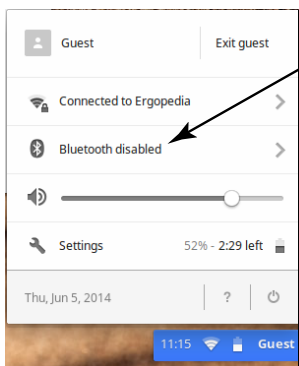


4. Turn Bluetooth on by clicking on "Enable Bluetooth."

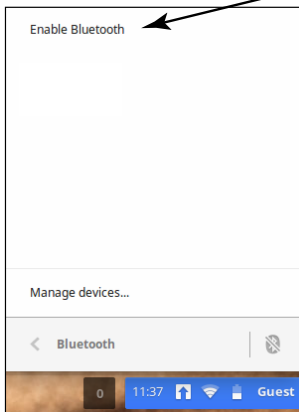

The computer will immediately look for available Bluetooth devices.

5. Select the ErgoBot device to which you want to pair and connect.

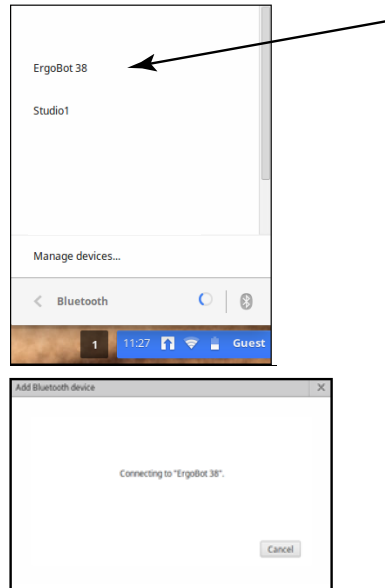

6. This shows a successful connection.



7. Sometimes you may have to turn the ErgoBot power switch on and off to refresh the Bluetooth connection.

You can now run the interactive elements in the *Essential Physics e-Book* with your paired ErgoBot!