


Pairing your computer to the ErgoBot: **Windows**

1. Turn on the ErgoBot using the switch on its bottom.

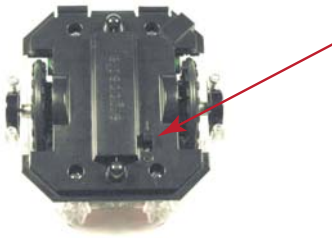

2. Click the **Bluetooth** button on the system tray in the lower right corner of your computer's Desktop.




3. Select **"Turn Adapter On"** to turn on Bluetooth if it is currently off. (You can also access this by clicking Windows -> Control Panel -> Devices and Printers.)

Add a Device
Show Bluetooth Devices
Open Settings
Turn Adapter On


4. Click the System Tray's Bluetooth button again and now select **"Show Bluetooth Devices."**

Add a Device
Show Bluetooth Devices
Open Settings
Turn Adapter On


5. Select **"Add a device"** to add the new ErgoBot as a Bluetooth device to your computer.

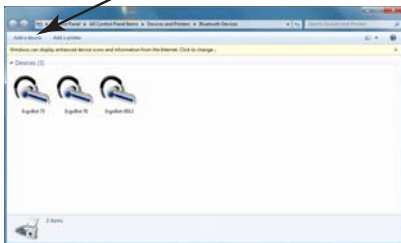

6. Select the device you wish to add and click **"Next."**

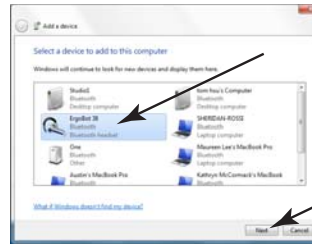


7. The device has now been added successfully to the computer. Click **"Close."**

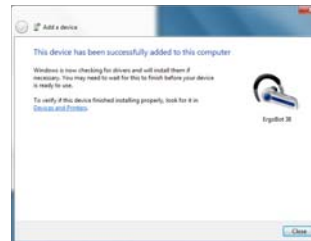

8. Double-click your new device in the control panel for Bluetooth devices.



9. This shows a successful connection.

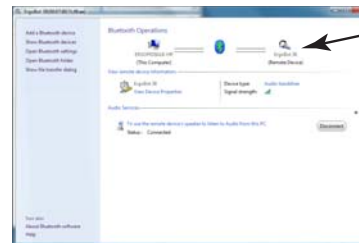

You can now run the interactive elements in the *Essential Physics* e-Book with your paired ErgoBot!